


FLOW CYTOMETRY

MoFlo™ 4Way™ Sort

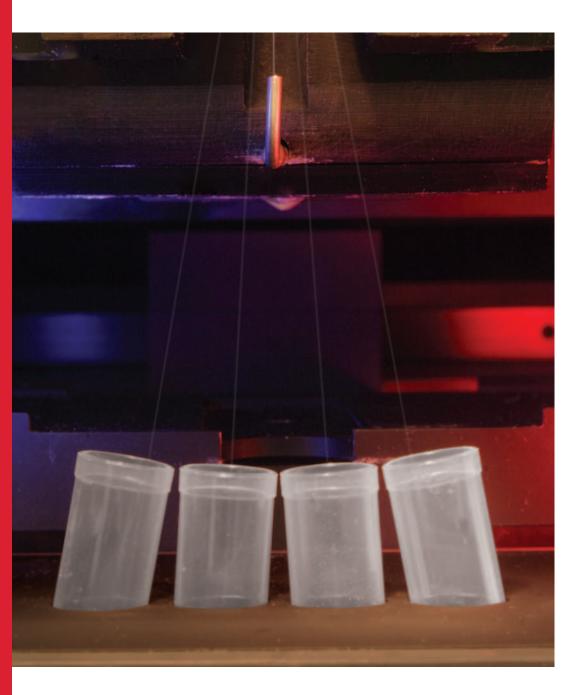

Power and flexibility for a fully integrated, automated sort.


The patented 4Way Sort system upgrades the standard two-stream capability of the MoFlo High-Performance Cell Sorter to a three-and four-stream capability.


Researchers can sort more subsets faster and with higher viability, maximizing the sample's purity and increasing productivity in your lab.


MoFlo 4Way Sort


Based on unlimited Boolean gate logic, the patented 4Way Sort system enables operators to define up to four unique populations of interest (Figure 1).

- Increase productivity and save precious sample volume with simultaneous sorting of four populations.
- Accommodate a variety of standard tube sizes and cuvettes.
- Manipulate charging for optimal droplet separation and maximum purity.
- Routinely switch between 2Way and 4Way sorting modes.
- 4Way Sort rates are not compromised by increased complexity of boolean gating logic.

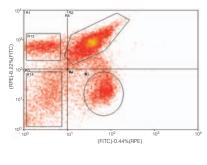

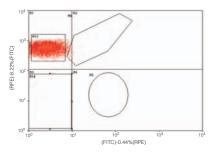

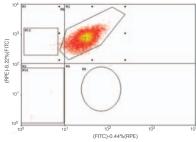

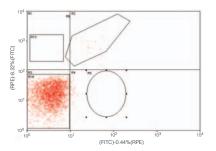

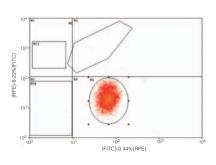


Figure 1. Mouse thymocytes were sorted into four subsets simultaneously with purity >99 percent in each tube.

4Way Sort Specifications

Collection Tube Sizes	1.5 mL, 5 mL, 7 mL
Stream Separations	13 mm
Maximum Cells per Second	70,000
Maximum Droplet Formation Rate	200,000 drops per second
Instrument Compatibility	SmartSampler, SortMaster™

For Research Use Only. Not for use in diagnostic procedures.
U.S. patent no. 5,483,469 Multiple Sort Flow Cytometer January 9, 1996.



Corporate Headquarters Denmark Tel +45 44 85 95 00

www.dako.com

Distributors in more than 50 countries

Australia +61 2 9316 4633

Austria +43 1 408 4334 50

Belgium

+32 016 38 72 20 Canada

+1 905 858 8510

Czech Republic +420 541 42 37 10

Denmark

+45 44 85 97 56 France +33 1 30 50 00 50

Germany +49 40 69 69 470 Ireland +353 91 768150

Italy +39 02 58 078 1

Japan +81 75 211 3655 The Netherlands +31 20 42 11 100 Norway +47 23 14 05 40

Poland +48 58-661 1879 Spain

. +34 93 499 05 06 Sweden +46 08 556 20 600 Switzerland +41 41 760 11 66

United Kingdom +44 (0)1 353 66 99 11

United States of America +1 805 566 6655