

SOP-P067

Rectangle Plate Procedure on back of sheet

Making Agar for Bacteria Growth Round Plates

Objective: To prepare Agar for growing bacteria for the Bardot Project

Procedure: Use the Trypticase Soy Agar Catalog #211043 BD (500g)

1. Weigh out 40gm of the powder and suspend it in to 1L of Millipore Water.
2. Mix thoroughly. Heat with frequent agitation and boil for 1 minute to completely dissolve the powder. Do not overheat. Then Autoclave.

Procedure: Pour Agar into Petri Dishes: From the refrigerator method.

1. Remove the bottles containing 500ml of the agar from the refrigerator and let it warm up to room temperature. Each bottle should fill 18 to 20 petri dishes.
2. Be sure you have a microwaveable container with at least 500ml of water in the microwave with the agar. Place the agar in the microwave with the lids loosened for 15 to 20 minutes on high. The agar must be completely dissolved so you may need to cook it longer.
3. Once dissolved, place the bottles in the pre-warmed water bath, set at 60°C. The agar should be around 55°C before pouring. This usually takes approximately 5 minutes.
4. Then remove from the water bath and place the bottles on the flat surface. Place the sterile petri dishes on the flat surface to be filled with the agar.
5. **Pipette 25ml** (draw up 27 but use only 25ml) of the agar into the petri dishes being careful not to create bubbles. If bubbles do occur, just remove them with the pipette. Let cool and then place them with the agar side up carefully in the plastic sleeve they came in. Seal the sleeve with tape and label each stack the type of agar and the date. The next day, place the plates you poured in the 37° incubator for the sterility test overnight. If no contamination, label each stack “sterility check-good”. Then store in the cold room until needed.

Procedure: Pouring plates of Agar without chilling first:

1. Once autoclaved, let the Agar cool down to safe handling temperature. (Bottles will still be warm to touch.) Place the bottles containing Agar in a 60°C waterbath.
2. Repeat steps 4 and 5 above.

Rectangle Plates Procedure on back ➔

Created by: Cheryl Holdman

Verified by: _____ **Date:** 10/22/2010 _____

Print Name

Sign Name