

SOP-M014

Nasty Bardot

Objective: To view Scatter Patterns of Various Bacteria Colonies

BE SURE YOU ARE ON THE CFSE DATABASE BEFORE YOU BEGIN.

1. Click on the Lady Bug Icon on the lower bar to open the Bug Buster Software.
2. Place the plate with Agar side down in the provided holder.
3. Click the “Take scan image” icon on the upper bar. The colonies will appear as red dots.




If there are any multiple groups or odd shaped colonies, you can use the “Hand” icon to turn the colonies from red to green. Now the Bardot will only scan the red ones.




If you want to start your scan with a particular colony, use the “Plus Sign” icon. Go to the specific colony you want and click once with the mouse.

4. Now you are ready to start scanning. Click on the “Take scatter pictures” icon.

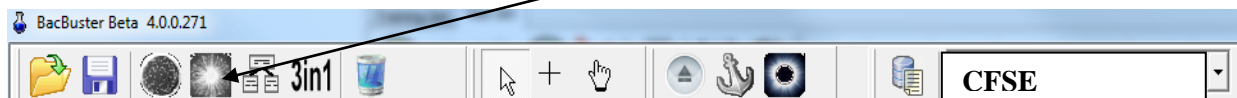

Saving Data:

Here is an example of the information needed when saving the data to the database.

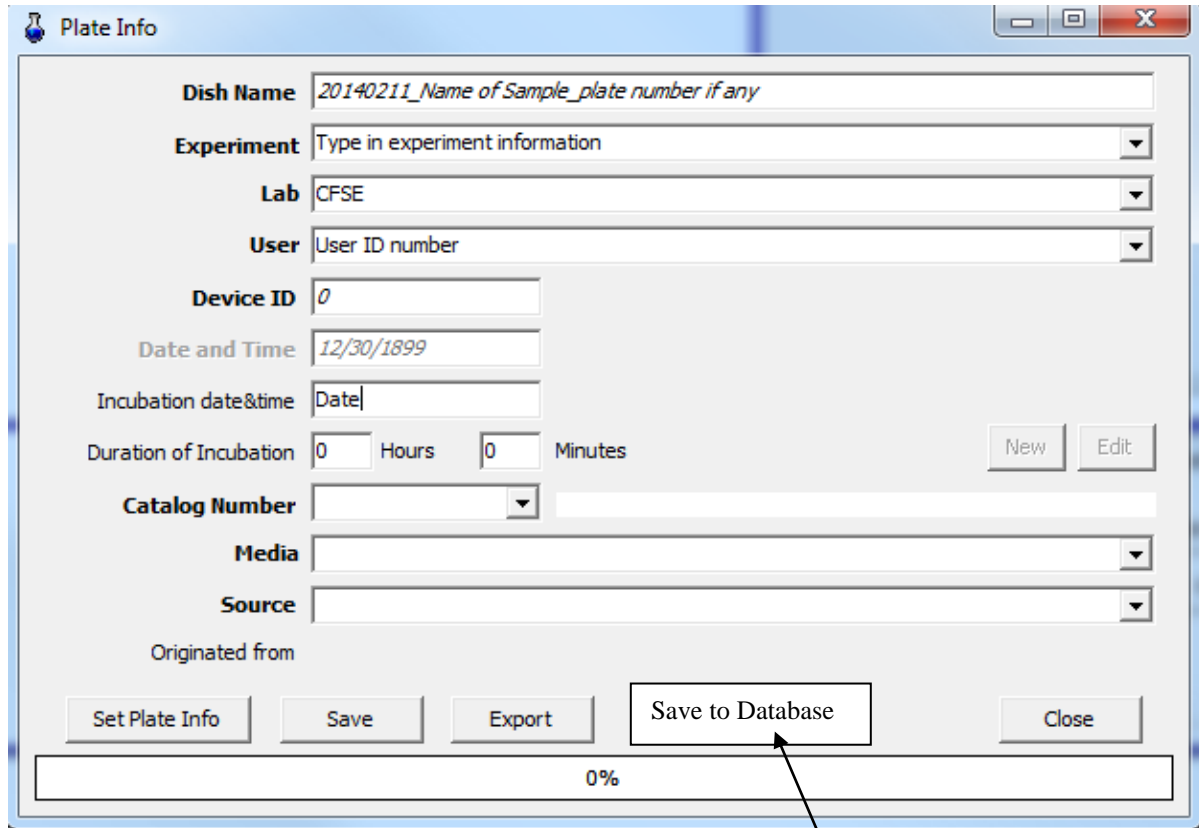


The screenshot shows a Windows-style dialog box titled "Plate Info". It contains several input fields and buttons. The fields are: "Dish Name" (text input with placeholder "20140211_Name of Sample_plate number if any"), "Experiment" (dropdown menu with "Type in experiment information"), "Lab" (dropdown menu with "CFSE"), "User" (dropdown menu with "User ID number"), "Device ID" (text input with "0"), "Date and Time" (text input with "12/30/1899"), "Incubation date&time" (text input with "Date"), "Duration of Incubation" (two spinners for "Hours" and "Minutes", both set to "0"), "Catalog Number" (dropdown menu), "Media" (dropdown menu), and "Source" (dropdown menu). There are also "New" and "Edit" buttons. At the bottom, there are buttons for "Set Plate Info", "Save", "Export", "Save to Database", and "Close". A progress bar at the very bottom shows "0%". An arrow points from the "Save to Database" button to the text below.

After all the information has been entered correctly, click the “Save to Database”


Note: The information MUST be entered in the same way each time.

Analysis:

Use the “Bacterium Analyzer5” software. Icon is on the lower information bar.

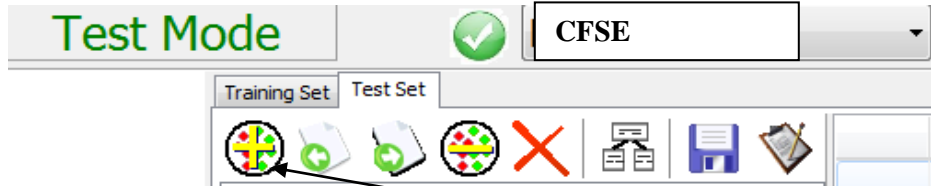

Be in Test Mode / Test Set Tab / and then click on the “Add Plate” icon. A window will open displaying your plate information.

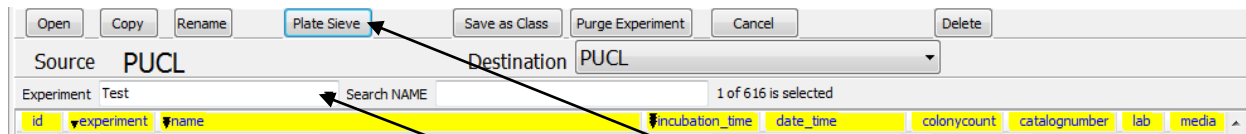

Using the drop down arrow, choose you experiment name.

If you have several plates and want to combine the data for comparison, use the “Plate Sieve” tab.

This could take a few minutes if there is a large amount of data.