## **SOP-P036**

## Making 3% Dextran

**Objective**: To make 3% Dextran to be used in the Dextran sedimentation procedure (SOP#-P035) for the isolation of neutrophils from whole blood.

## **Procedure**:

- 1. Measure out 3gm of Dextran T500 (Pharmacia #17-0320-01).
- 2. Combine in 150ml flask with 100ml of Phosphate Buffered Saline(PBS). See SOP#-P020 for preparation of PBS.
- 3. When Dextran is dissolved (no particulate is seen) transfer to a proper storage container, such as a 100ml Wheaton bottle with lid.
- 4. Label the bottle with contents, date made, and initials.
- 5. Store at 4°C.

Created by: Kathy Ragheb	DATE: <u>2/25/99</u>	
Verified by:	Date:	
Print Name		Sign Name